



emulate

EP-244 Chip-Array Basic Research Kit Protocol

Rev B

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 2 OF 64

Experimental Overview

- **Equipment and Materials Required**
- **Workstation Preparation, Chip-Array Handling Techniques**
- **Pre-Experiment: Reagent Preparation**
- **Prior to Day -2: AVA Setup**
- **Day -1: Chip-Array Preparation**
 - Prepare Chip-Arrays
 - Prepare ECM solution
 - Coat chips with ECM
- **Day 0: Introduce Cells to Chip-Arrays**
 - Prepare Chip-Arrays
 - Prepare Cells
 - Seed cells to top channel
- **Day 1: Chip-Array Prime, Regulate, and Flow**
 - Wash Chip-Arrays
 - Seed cells to bottom channel
 - Gas equilibration of media
 - Prime Chip-Arrays
 - First Regulate Chip-Arrays, Flow
- **Day 2: Chip-Array Regulate**
 - Second Regulate Chip-Array, Flow
- **Day 5 +: Chip-Array Maintenance and Sampling**
 - Sampling and media replenishment
 - Routine automated imaging of biology
 - Dosing
- **Troubleshooting**

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 3 OF 64

Equipment and Materials Required

Introduction

Ensure all equipment, materials, and reagents are accessible prior to each experiment. Below are the catalog numbers for the specific materials needed.

Note: Exact catalog numbers are not provided for some required materials, as several brands and models are accepted. For those materials with specific catalog numbers we recommend purchasing this exact item in order to guarantee success.

Required Emulate Products

A list of Emulate products needed for this protocol is provided below:

Item	Quantity/Description	Vendor	Cat #
AVA™ Emulation System	1 Instrument	Emulate	AVA-ES1
Orb-HM1® Hub Module	1 Module	Emulate	ORB-HM1
Chip-Array Basic Research Kit-96-Emulation Pack	• Chip-Array Organ-Chip consumables (8) • Pipette Guides (3 per Chip-Array; 24 total) • Flow Guides (3 per Chip-Array; 24 total) • Pipette Guide Covers (1 per Chip-Array; 8 total) • Steriflip (14 total)	Emulate	BRK-CA-96

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 4 OF 64

Required Lab Equipment and Consumables

A list of lab equipment and consumables needed for this protocol is provided below:

Item	Quantity/Description	Vendor	Cat #
MycoFog® Biodecontamination Starter Kit	1	MycoFog®	MF-SK-1Bx
Steriflip-HV Filters (Sterile, 0.45 µm PVDF-filter)	25/pack	EMD Millipore	SE1M003M00
Handheld vacuum aspirator	1 or 2 aspirators for BSC	Corning	4930
Aspirating pipettes (2-mL, polystyrene, individually wrapped)	-	VWR	53106-450
Aspirating tips	-	-	-
Serological pipettes (2-mL, low- endotoxin, sterile)	-	VWR	55300-363
Serological pipettes (5-mL, low- endotoxin, sterile)	-	VWR	55300-421
Serological pipettes (10-mL, low- endotoxin, sterile)	-	VWR	55300-523
Serological pipettes (25-mL, low- endotoxin, sterile)	-	VWR	55300-567
Serological pipettes (50-mL, low- endotoxin, sterile)	-	VWR	53106-441
P20 Single Channel Pipette	-	-	-
P200 Single Channel Pipette	-	-	-
P1000 Single Channel Pipette	-	-	-

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 5 OF 64

P20 Multi Channel Pipette	-	-	-
P200 Multi Channel Pipette	-	-	-
P1000 Multi Channel Pipette	-	-	-
P20 sterile, low-adhesion, filtered pipette tips*	-	VWR	76322-134
P200 sterile, low-adhesion, filtered pipette tips	-	VWR	76322-150
P1000 sterile, low-adhesion, filtered pipette tips*	-	VWR	76322-154
15 mL Falcon Conical Tubes	-	VWR	21008-918
50 mL Falcon Conical Tubes	-	VWR	21008-940
50 mL Disposable Pipetting Reservoirs	5/pack	VWR	89094-682
Nunc™ 96-Well Polystyrene Round Bottom Microwell Plates	10/pack	ThermoFisher	268200
Nunc™ 96-Well Polypropylene DeepWell™ Sample Plates, Square, Conical Bottom	50/pack	ThermoFisher	95040462
Eppendorf tubes	-	-	-
Aluminum foil	-	-	-
Parafilm M Sealing Film	-	VWR	52858-076
Microscope (with camera), for phase contrast imaging	-	-	-
Hemocytometer	-	-	-

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 6 OF 64

Water bath (or bead bath)	-	-	-
Vacuum set-up	-	-	-
Centrifuge	-	-	-
Ice bucket	-	-	-
70% ethanol	-	-	-

***Disclaimer: The listed P20 and P200 filtered pipette tips are recommended due to their validated robustness in interfacing with the Chip-Array.**

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 7 OF 64

Workstation Preparation and Chip-Array Handling Techniques

Overview

Aseptic Techniques

- Always work with chips in a sterile environment, such as the biosafety cabinet (BSC).
- Before beginning the experiment, prepare, sanitize, and organize the work surface of the BSC. Arrange tips, media, and other necessary materials in the sterile field so they are easily within reach but not blocking the path of airflow.
- Use 70% ethanol to thoroughly sanitize all reagents and materials before placing them into the BSC.
- Always avoid touching the Chip-Array Emulations directly. Handle the Chip-Arrays only by the sides, by the lid, or by the tab, with gloves.

Cell Storage

Always store cryopreserved cells in the vapor phase of liquid nitrogen. Never store them in dry ice or in a -80°C freezer. Chronic temperature fluctuations can cause severe damage to cell membranes and cytoskeletal components.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 8 OF 64

Chip-Array Handling Techniques

Chip-Array Design

Figure 1 shows the Chip-Array. It can be divided into two separate sections: the “Reservoir” section and the “Emulations” section.

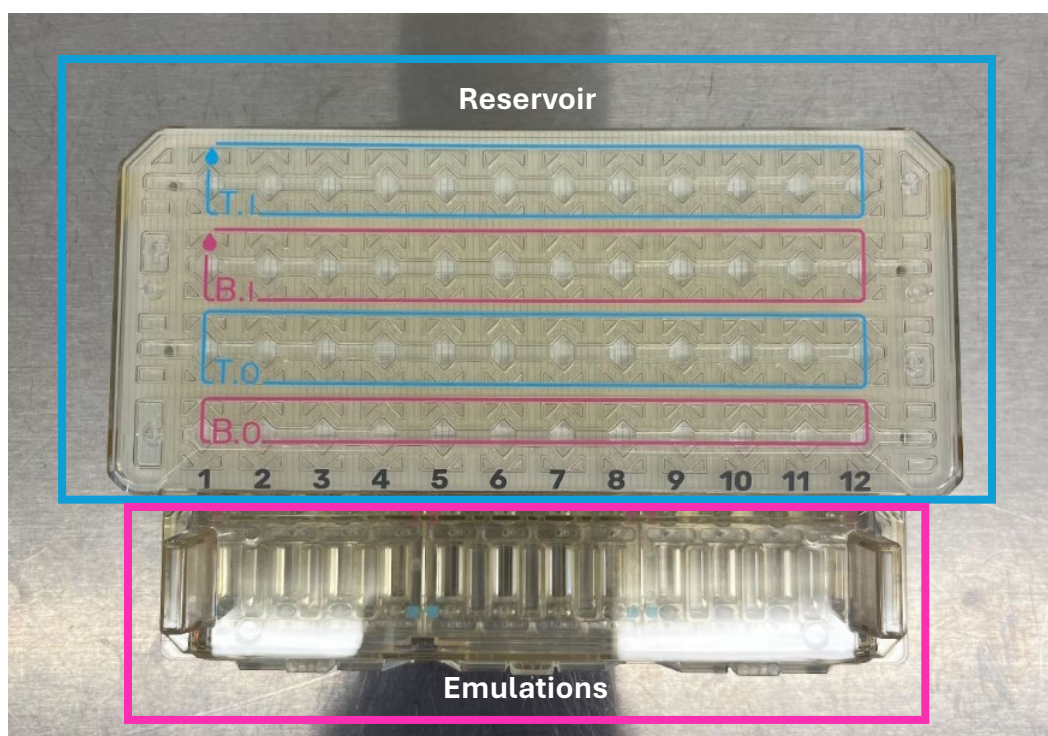


Figure 1. Chip-Array consumable design.

One Chip-Array contains twelve “Emulations” (see Figure 2 below). An Emulation is an independent Organ-Chip sample; each Emulation has its own top and bottom microfluidic channel, which is separated by a porous membrane. Each Emulation is compartmentalized from one another, meaning that each Emulation can be treated individually from each other. Each Emulation is labeled 1-12 on the lid of the reservoir.

Additionally, each Emulation has its own set of 4 reservoirs: a top channel inlet reservoir (TI), a top channel outlet reservoir (TO), a bottom channel inlet reservoir (BI), and a bottom channel outlet reservoir (BO). This is indicated by the label on the reservoir lid.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 9 OF 64

The inlet reservoirs are where media is added during media replenishment and dosing, which will be supplied to the biology seeded in the Emulations via unidirectional flow facilitated by the AVA instrument. Because each Emulation's reservoirs are compartmentalized from each other, different compounds and concentrations can be dosed into each Emulation independently from one another, allowing users to have multiple dosing conditions on a single Chip-Array.

Each Emulation can be broken down into separate inlet and outlet ports for the top and bottom channels. The inlet sides of each channel are towards the reservoir, while the outlet sides are oriented away from the reservoir. Inlet and outlet orientation are based on the direction of flow once the Chip-Array is under flow in the AVA instrument. Media will flow unidirectionally from inlet to outlet, from the inlet reservoir through the Emulation then into the outlet reservoir.

Additionally, a viewing window is cut into the gasket on each Emulation, allowing users to image directly into the channels. Each Emulation is labeled with its own chip number (1 through 12) and is visible under the microscope.

Chip-Arrays come with tape on the reservoir lid. This helps to prevent the lid from accidentally falling off when users are handling the Chip-Array and any inversion steps when seeding the bottom channel. This tape can be removed once users are ready to start adding media to the reservoir during the Prime step.

See Figure 2 for a graphic that displays the orientation of each Emulation.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 10 OF 64

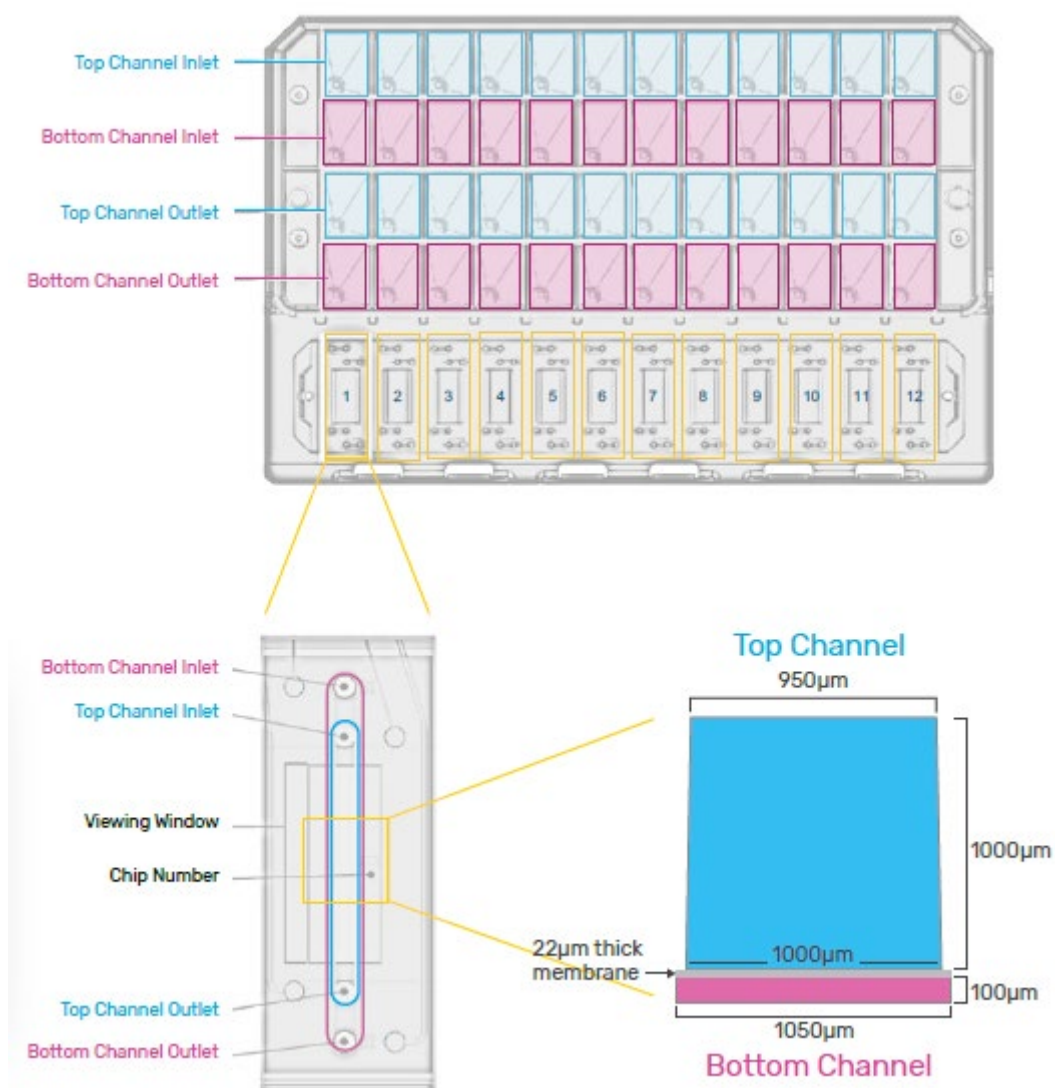


Figure 2. Schematic of an Emulation within the Chip-Array consumable.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 11 OF 64

Components of the Chip-Array

One Chip-Array will come in its package with three pipette guides, three flow guides, and one pipette guide cover. See Figure 3 for what these Chip-Array components look like.

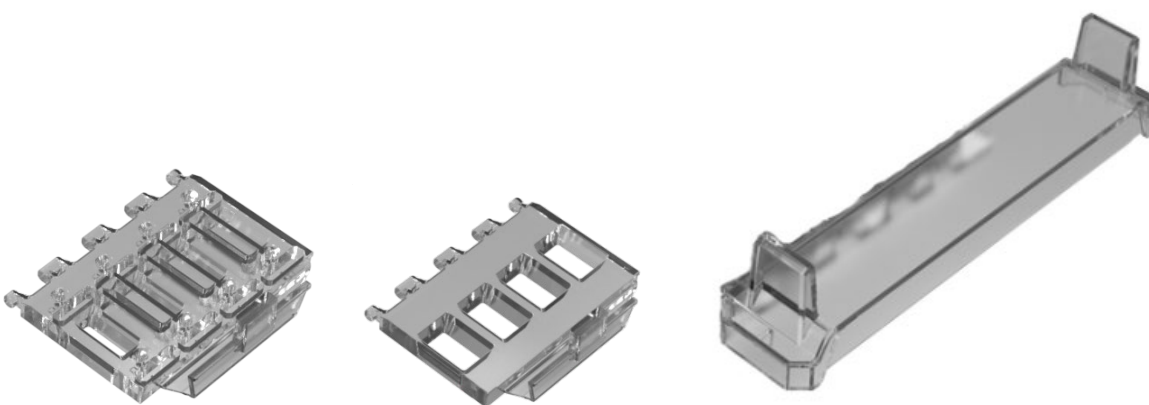


Figure 3. Chip-Array accessories: pipette guide (left), flow guide (middle), and pipette guide cover (right).

Pipette Guides

Pipette guides facilitate manual pipetting with micropipettes into the top and bottom channels of each Emulation on the Chip-Arrays. These channels are separated by a porous membrane and therefore have separate pipetting/seeding ports that allow users to pipette into each channel individually. Single channel and multichannel pipettes are compatible with the Chip-Array pipette guides. Each pipette guide interfaces with 4 Chip-Array Emulations, meaning that one Chip-Array accommodates three pipette guides.

For each Emulation, the pipette guide contains two seeding ports, which allow the user to manually pipette solution and cell suspension into the top and bottom channels using a single channel or multichannel pipette. **These seeding ports are compatible with P20 and P200 pipette tips. Refer to the Required Lab Equipment and Consumables Section for recommended pipette tips.** The top channel seeding port is at the bottom of the pipette guide, while the bottom channel seeding port is at the top of the pipette guide. On each pipette guide, the magenta color indicates the location of

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 12 OF 64

the seeding ports used for pipetting into the bottom channel, while the blue color indicates the location of the seeding ports used for pipetting into the top channel.

Additionally, the pipette guide contains outflow troughs for each channel, which collects the outflow from each channel as users pipette into them. The bottom channel trough is on the opposite side of the bottom channel seeding port, while the top channel trough is on the opposite side of the top channel seeding port. These troughs can be filled with excess cell culture media or extracellular matrix (ECM) solution when pipetting in order to prevent evaporation of the liquid in the Emulations during long periods of incubation (e.g., overnight).

Pipette guides also contain humidity PBS reservoirs between each Emulation (3 humidity reservoirs per pipette guide). These reservoirs should be filled with PBS to ensure minimal evaporation of liquid in the Emulation channels.

See Figure 4 for the orientation of the different ports and wells of the pipette guide.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 13 OF 64

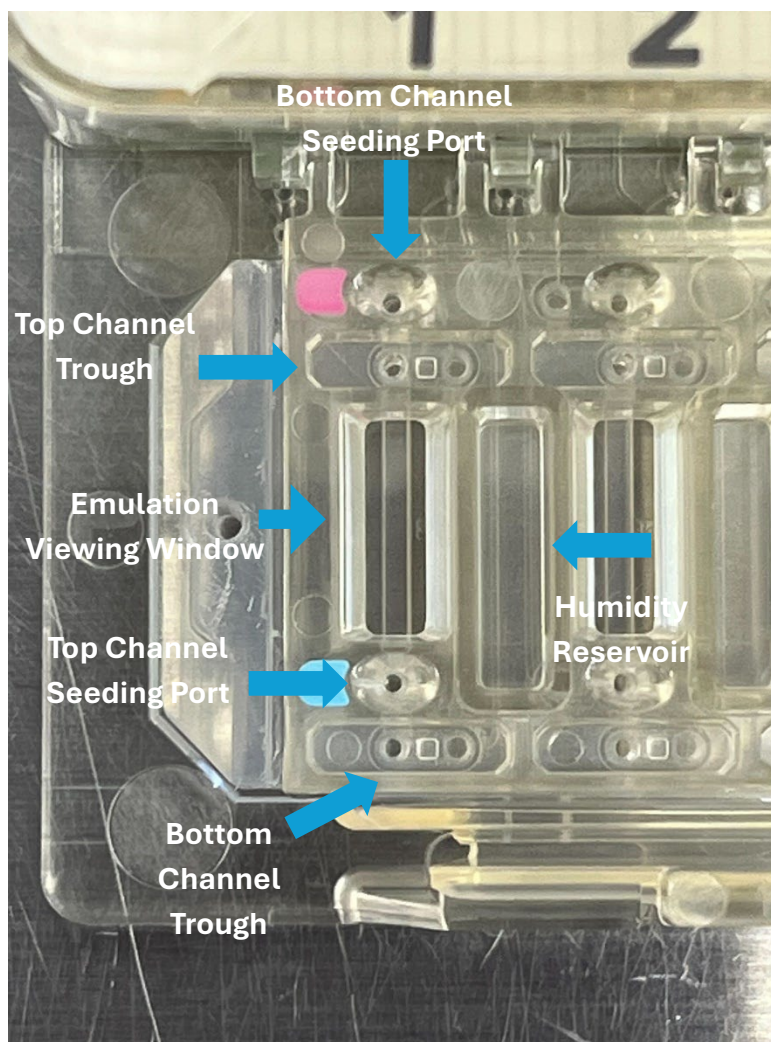


Figure 4. Orientation and components of the pipette guide.

Pipette Guide Covers

Pipette guide covers act as a shield to help maintain sterility while pipette guides are attached to the Emulations. Pipette guide covers should always be used when pipette guides are attached to the Chip-Array and should only be removed when users are handling the Chip-Array in the BSC. The covers are not required once flow guides are attached to the Chip-Array, as the flow guides provide a complete seal that prevents exposure of biology in the culture channels to the outside environment.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 14 OF 64

It should be noted that pipette guide covers have a specific orientation. The larger, flat side of the pipette guide cover should face the user, while the shorter, angled side should face the Chip-Array reservoir. This specific orientation allows the pipette guide cover to properly slot over the Emulations and remain flush. See Figure 5 for the proper orientation of the pipette guide cover.

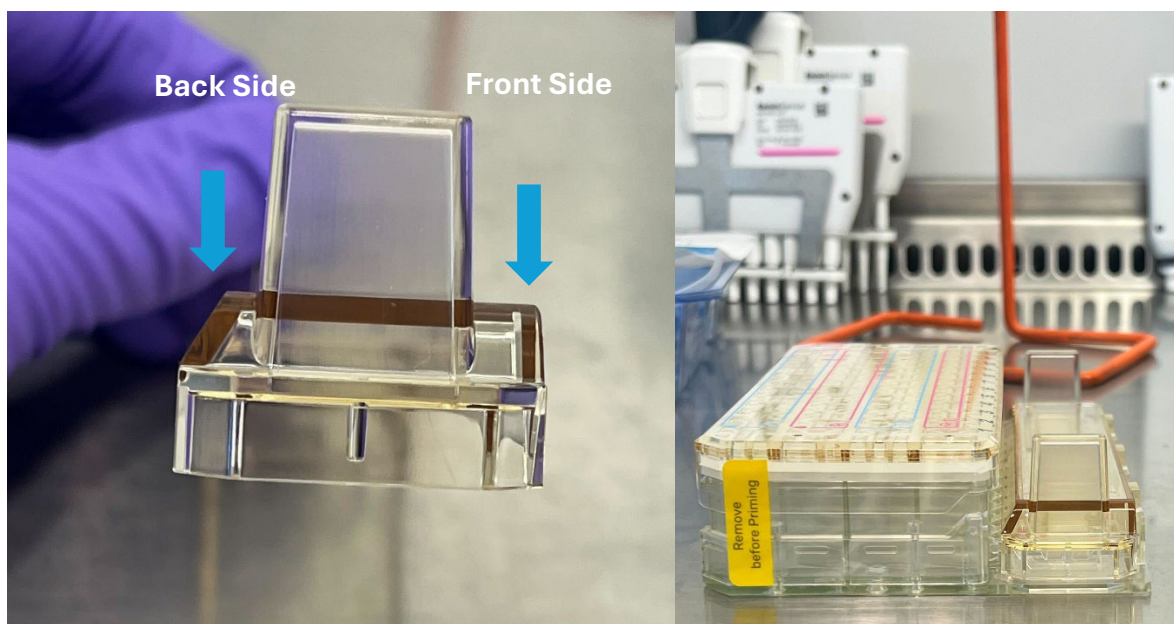


Figure 5. Proper orientation of the pipette guide cover.

Flow Guides

Flow guides are used to fluidically connect the Emulations to the reservoirs filled with media, allowing the AVA to flow media from the reservoir inlets to the Emulations.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 15 OF 64

Pipetting Solution into Channels

All wash steps are typically performed with a P200 multichannel pipette, using 50-100 μ L of the specific solution, which is enough to fill the channels and leave excess in the troughs. While the Chip-Array is compatible with multichannel pipettes using 12 tips at a time, it is recommended for users to use 6 tips at a time during these wash steps. When using multiple tips at a time, users must ensure that all tips are properly interfacing with the seeding port, and enough force is applied to the tips. If there is not a proper seal, liquid will not flow through the channel.

Follow the steps below for pipetting solution into each Chip-Array using a multichannel P200 pipette.

1. Take a multichannel P200 pipette with up to 6 sterile pipette tips and collect the solution to be added to the Chip-Array Emulations. Ensure that all tips on the pipette are properly secure and attached. Failure to do so can cause tips to fall off during handling or for tips to collect the incorrect volume of solution.
2. Align all pipette tips with the specific channel seeding port and insert all tips perpendicular to the port, ensuring that the tips are secure in the port.
 - a. Users must ensure that all tips are properly secured in the seeding port and enough force is applied by the user during insertion of tips. Lack of a proper seal will cause improper liquid dispensing.
 - b. When pipetting with multiple tips, users can use their offhand to steady the pipette with their fingers on the pipette shoulders to distribute force before pipetting. Always make sure tips are perpendicular when inserted into the pipette guide; failure to do so can cause tips to become stuck.
3. Steadily dispense the liquid through the channel. See Figure 6 for proper placement of tips into seeding port and dispensing of liquid through the channels.
4. Excess liquid will flow into the trough of the channel that is being washed. Users can either keep the excess liquid in the through or aspirate the liquid.
5. If aspirating liquid from the troughs, users must avoid touching the port in the trough, as this will aspirate the liquid that is in the Emulation channels. It is recommended for users to place their aspirating tips along the left or rightmost walls of the trough that is being aspirated. See Figure 7 for proper placement of tips when aspirating from troughs.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 16 OF 64

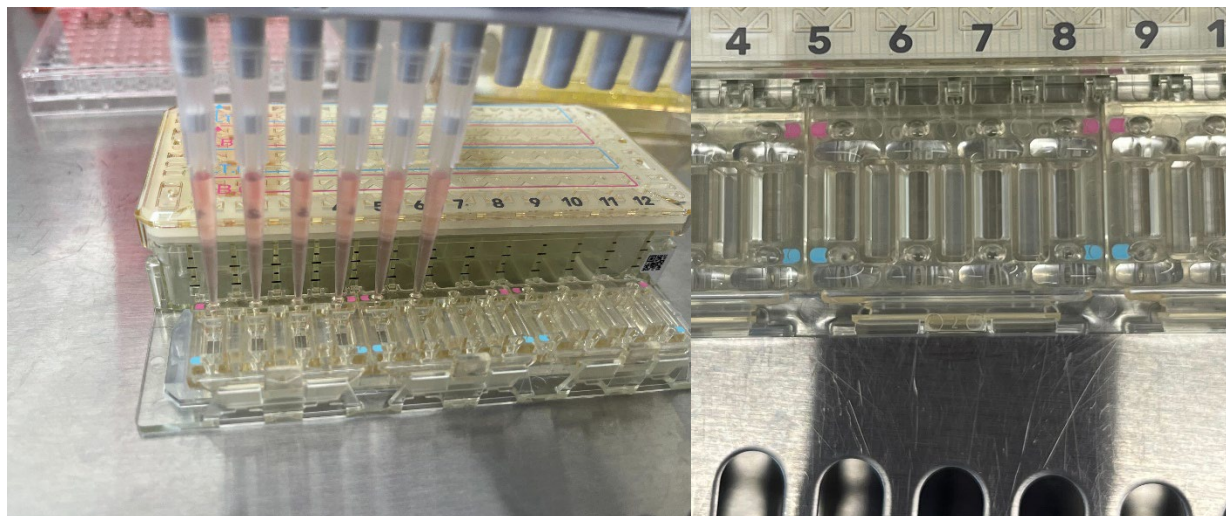


Figure 6. Left image shows proper placement of tips into seeding ports and dispensing of liquid. Right image shows collection of excess outflow into the trough.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 17 OF 64

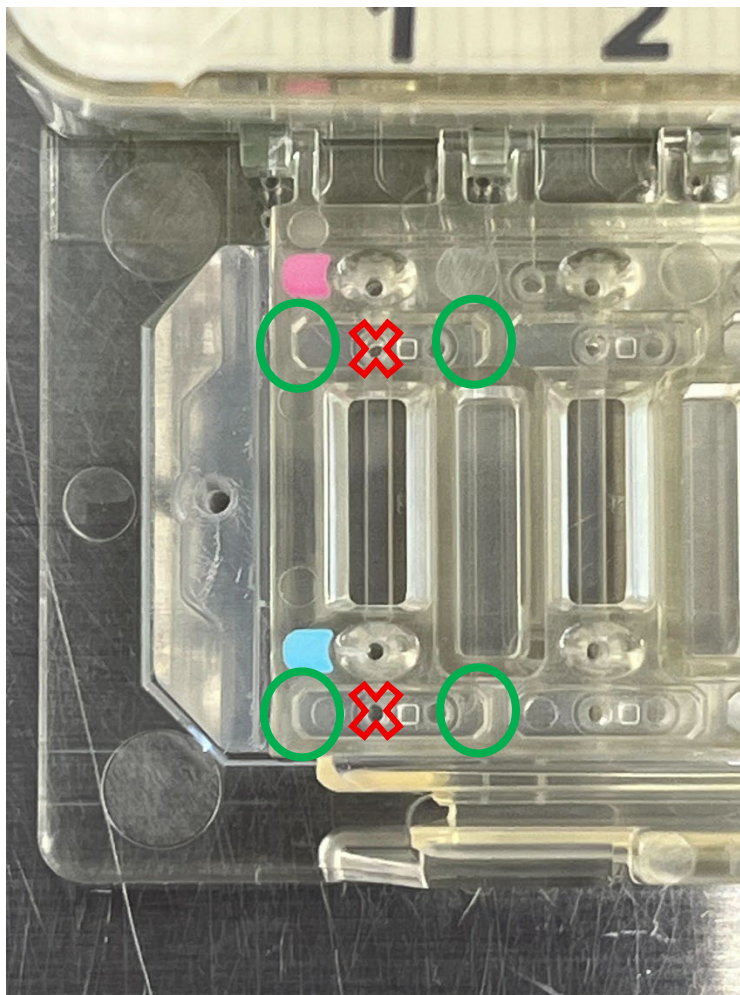


Figure 7. Proper placement of tip when aspirating from trough.

CAUTION: If excess liquid during pipetting falls into the viewing window of the Emulation, take precaution NOT to make direct contact with the viewing window with the aspirating pipette tip. The top gas exchanger film can be easily punctured by a pipette tip. Tips should be placed in the corner of the viewing window and gently lowered until all excess liquid is aspirated from the viewing window.

All seeding steps are performed with a P20 pipette using the same steps above. **It is recommended that users seed a maximum of 6 Emulations at a time. Seeding with more than 6 at a time increases the likelihood of variability between**

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 18 OF 64

Emulations. The table below highlights the seeding volumes recommended per channel.

Channel	Volume
Top Channel	20 μ L
Bottom Channel	10 μ L

CAUTION: Users must be careful when using P20 pipette tips during cell seeding steps. P20 pipette tips are more likely to get stuck in the seeding ports during seeding compared to P200 tips during washing. If a tip gets stuck, users can pull the tip out of the seeding port. However, it is likely that this will compromise cell distribution within the channel of that specific Emulation, likely leading to a re-seed of that Emulation. Users should follow these guidelines to prevent tips getting stuck:

- Always make sure the pipette tips and multichannel pipette are perpendicular to the seeding port and vertical. Removing tips from the seeding ports at an angle can increase likelihood of the tips getting stuck.
- When seeding several Emulations at a time with P20 pipette tips, users must be careful with the amount of force applied to the multichannel P20 tips. Inserting the tips is enough to create a seal with the seeding port to ensure proper flow. Additional excessive force when pipetting multiple Emulations at a time (i.e. using your offhand to apply force to the multichannel pipette) can cause tips to become stuck in the pipette guide seeding port.
- Refer to the Required Lab Equipment and Consumables Section for recommended pipette tips. These pipette tips are recommended as the tip shape is most compatible with the pipette guide seeding ports and has been thoroughly tested.

CAUTION: When seeding, take care to aspirate all excess liquid from the trough of the specific channel that is being seeded. This should be done before and after pipetting cells through the specific channel seeding port. Any excess liquid in the trough may contribute to improper cell distribution and gradients during seeding.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 19 OF 64

NOTE: When seeding, it is recommended to leave the P20 pipette tip in the seeding port until the outflow volume in the trough is aspirated, then the user may remove the pipette tip.

Attachment and Detachment of Pipette Guides and Flow Guides onto Chip-Array

There are three sections of the Chip-Array that allow connection with a pipette guide or flow guide: left, right, and center. The left pipette guide or flow guide will interface with Emulations 1-4, the center pipette guide or flow guide with Emulations 5-8, and the right pipette guide or flow guide with Emulations 9-12. This means that each Chip-Array accommodates three pipette guides or flow guides at a time. See Figure 8 for a breakdown of these sections.

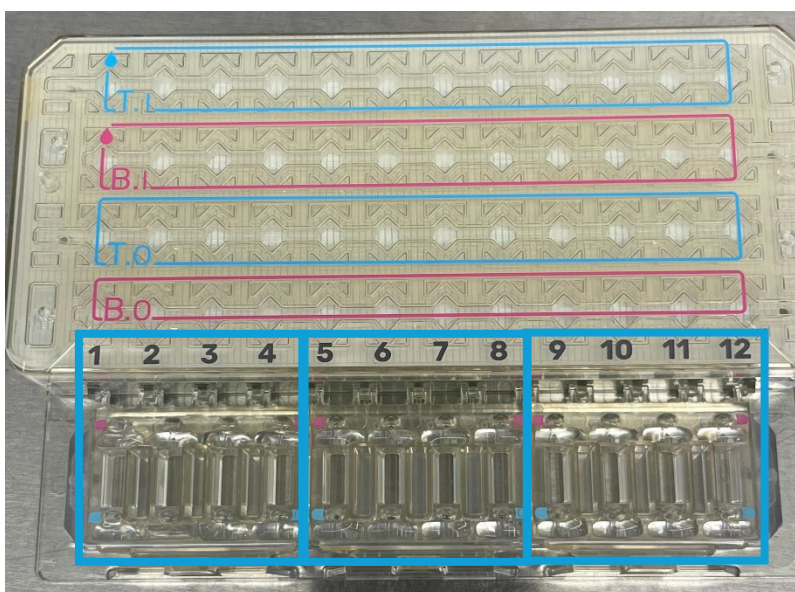


Figure 8. The three different sections of the Chip-Array for pipette guides or flow guides.

All Chip-Arrays will come packaged with three pipette guides attached to the consumable, meaning users will not have to attach the pipette guides and can immediately begin pipetting solution into the Emulations. Pipette guides should be detached and re-attached if any leakages are observed during ECM coating/washing steps. The following steps, while particular for flow guide attachment, are also applicable to pipette guide attachment.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 20 OF 64

Each guide is connected to the Chip-Array via hinges on the Chip-Array. The guides have connection points that interface with these hinges. See Figure 9, which highlights the connection points on the guides and the hinges on the Chip-Array.

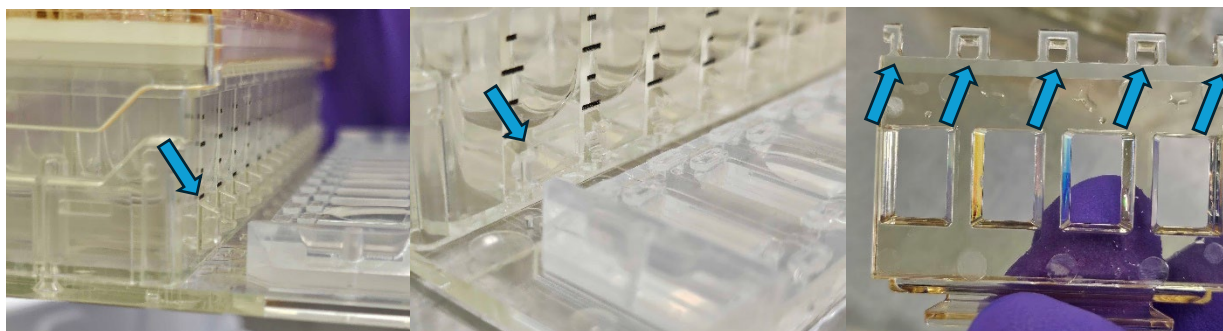


Figure 9. Pipette or flow guide connection points and hinges on Chip-Array.

It is recommended to attach the left guide first, followed by the middle guide, then lastly the right guide. Follow the steps below:

1. Align each guide connection point with the hinges on the Chip-Array. Push these connection points into the hinges; the user should be able to feel or hear a very light snapping noise, which indicates proper connection.
2. Once connected, gently snap the flow guides in place one at a time, using two fingers to hold onto the flow guide tab while pushing downwards, so that it interfaces with its respective Emulations. Users should hear an audible click when snapping the guide into place over its Emulations. See Figures 10 and 11 for visuals on connected guides and then snapping them into place
3. For the left guide, make sure to align the guide as flush as possible to the leftmost hinge connector. When connecting the right guide, make sure to align the guide as flush to the rightmost hinge connector as possible. When connecting the middle guide, use the middle hinge on the guide and the middle hinge connector on the Chip-Array as reference points.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 21 OF 64

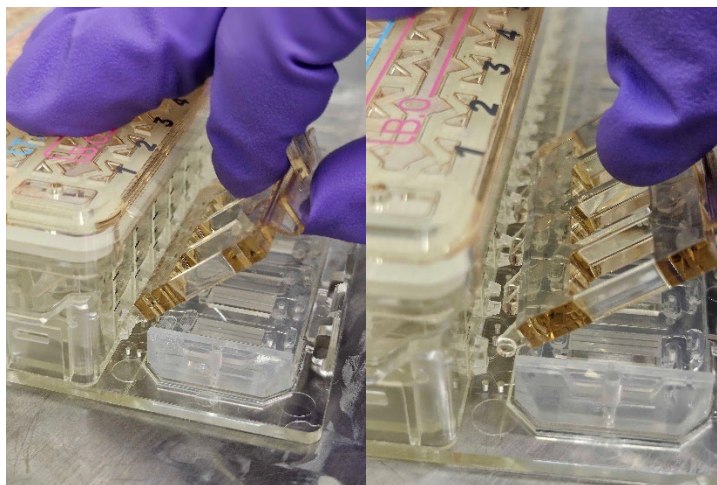


Figure 10. Connecting pipette or flow guide onto Chip-Array.

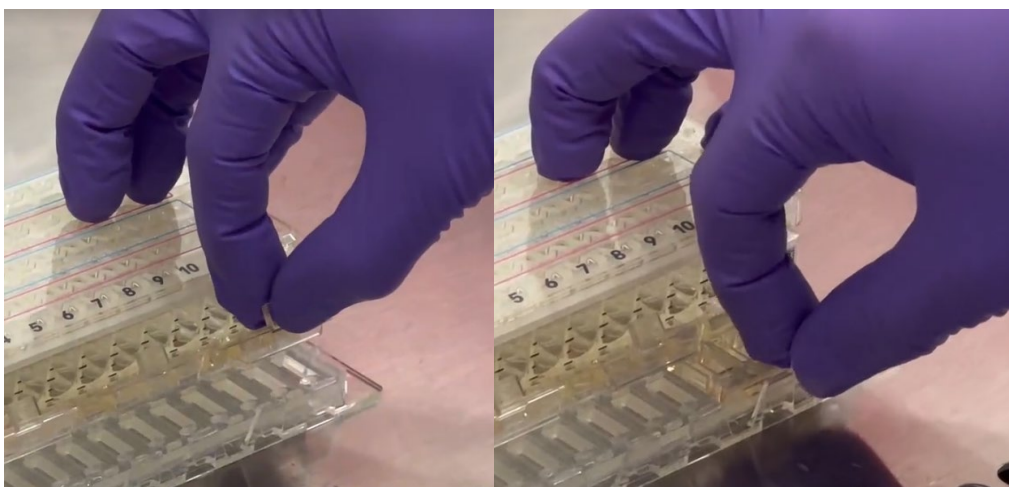


Figure 11. Snapping pipette or flow guide into place over Emulations.

Pipette and flow guides can also be removed from the Chip-Arrays after attachment. All Chip-Arrays come packaged with pipette guides attached to the consumables. Pipette guides should only be removed either if a leak is observed during pipetting, as this may indicate misalignment of the pipette guide with the Emulation, or during the Prime step which requires removal of pipette guides for the attachment of flow guides.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 22 OF 64

Remove guides by pushing on the tab that secures the guides inwards, then lifting the guide upwards. Once the clip is at a 45-degree angle, lightly tug on the clip to remove it. See Figure 12 for removal of the pipette guide.

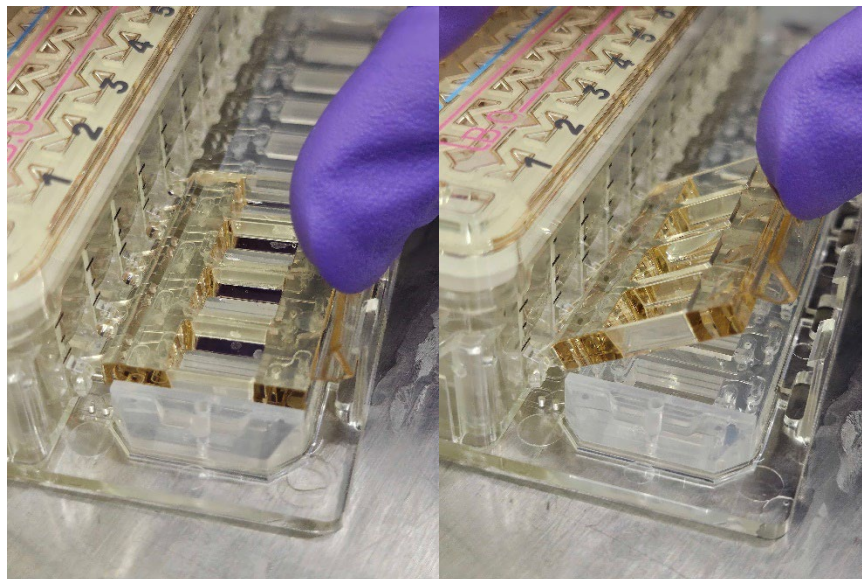


Figure 12. Removal of the pipette guide.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 23 OF 64

Prior to Day -2: AVA Setup

Overview

Goals

- Prepare the AVA instrument and software
-

AVA Decontamination

It is recommended by Day -2 to have the AVA decontaminated via MycoFog® after a service visit, or after a suspected contamination. This is approximately a 4-4.5 hour activity that requires heating the AVA to 37°C, then decontaminating via MycoFog®. If necessary, please contact Emulate Support for additional information about this process.

AVA Environmental Setup

After decontamination, it is recommended to have the AVA set to these environmental parameters:

- Temperature: 37°C
- CO2: 5%
- Ensure the water pan is filled to the 100% fill volume

Ensure that the environmental conditions of the AVA reach these specific setpoints and remain stable by Day 0.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 24 OF 64

AVA Experimental Design Setup

Steps

1. Open the Experiment Tab and create a new experiment using the Blank Template
2. In the Conditions window, input Experiment Name, Start Date of study and ensure that all cell, media, reagent, and ECM lot information is updated to reflect the materials used for the study
3. Input any necessary conditions, along with relevant information and number of replicates associated with each condition
4. In the Protocol window, input the desired protocol reminders and automated imaging routines for desired experimental days.
 - a. For example, schedule an automated imaging routine to collect images for cell health assessment. You can modify or add additional scheduled imaging routines at your discretion
5. Input the number of Chip-Arrays required, assigning names and serial numbers to each Chip-Array
6. Ensure experiment days in calendar view correctly align with expected timeline of activities

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 25 OF 64

Day -1: Chip-Array Preparation

Overview

Goals

- Coat the inner channels of Chip-Arrays with ECM proteins for cell attachment
-

Required Materials

- Chip-Arrays
 - 15 mL conical tubes
 - DPBS (-/-) at 4°C
 - ECM A
 - ECM B
 - 50 mL disposable pipetting reservoirs
 - 70% ethanol
 - Ice and ice bucket
 - Single channel and multichannel pipettes and filtered tips
 - Aspirator and sterile aspirating tips
 - Parafilm
-

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 26 OF 64

Chip-Array Preparation

Steps

1. Spray the Chip-Array packaging with 70% ethanol and bring into the BSC.
2. Open the packaging and take out all contents. This should include one Chip-Array, three pipette guides, three flow guides, and one pipette guide cover.
 - a. Pipette guides facilitate manual pipetting with micropipettes into the top and bottom channels of each Emulation on the Chip-Arrays
 - i. Single channel and multichannel pipettes are compatible with the Chip-Array pipette guides.
 - b. Flow guides are used to fluidically connect the Emulations to the reservoirs filled with media, allowing the AVA to flow media from the reservoir inlets to the Emulations.
 - c. Pipette guide covers act as a cover to help maintain sterility while pipette guides are attached to the Emulations. These are not necessary once flow guides are attached, as the flow guides provide a complete seal that prevents exposure of biology in the culture channels to the outside environment.
3. Chip-Arrays will already come with three pipette guides attached, so there is no need to attach the pipette guides.
4. Label each Chip-Array with its corresponding ID or label.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 27 OF 64

ECM Preparation

Before you begin

Prepare fresh ECM solution before each use by combining the individual ECM components (referred to as ECM A and ECM B in this protocol) with cold DPBS to reach the final working concentrations. The ECM solution will be used to coat both top and bottom channels. Prepare and maintain on ice.

Needed volumes

For human Chip-Arrays, the ECM working concentrations are:

Reagent	Final Concentration
ECM A	25 µg/mL
ECM B	100 µg/mL

Steps

1. Bring a full ice bucket to the BSC.
2. Calculate the volume of ECM solution needed to coat all Emulations.
 - a. Volume required per Emulation: ~140 µL
 - b. For every Chip-Array, prepare 2 mL of ECM solution:
 - i. 12 Emulations x 140 µL/Emulation = 1.68 mL of ECM solution
 - ii. 1.68 mL + extra 320 µL = 2.0 mL of ECM solution
3. Thaw the necessary number of aliquots of ECM A and ECM B on ice. Always maintain each ECM component and mixture on ice.
4. Combine the components to prepare the ECM working solution. See Example ECM Calculation section on next page for example of calculating ECM volumes.
5. Keep the ECM solution on ice until it is used.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 28 OF 64

Example ECM Calculation

ECM Calculation example for 8 Chip-Arrays

Solution	Concentration
ECM A stock concentration	1 mg/mL (C1)
ECM A final concentration	0.025 mg/mL (C2)
ECM B stock concentration	11.61 mg/mL (C1)
ECM B final concentration	0.1 mg/mL (C2)
Stock Volume	ECM A (X) or ECM B (Y) (V1)
Final volume of ECM solution	16 mL (V2)

ECM A Calculation:

$$(1 \text{ mg/mL}) \times (X \text{ mL}) = (0.025 \text{ mg/mL}) (16 \text{ mL})$$

$$Y = 400 \text{ } \mu\text{L of ECM A stock solution}$$

ECM B Calculation:

$$(11.61 \text{ mg/mL}) \times (X \text{ mL}) = (0.1 \text{ mg/mL}) (16 \text{ mL})$$

$$X = 137.81 \text{ } \mu\text{L of ECM B stock solution}$$

DPBS Calculation:

Volume DPBS =

$$(\text{total volume of ECM needed}) - (\text{volume of ECM A}) - (\text{volume of ECM B})$$

$$= 16000 \text{ } \mu\text{L} - 400 \text{ } \mu\text{L} - 137.81 \text{ } \mu\text{L}$$

$$= 15462.19 \text{ } \mu\text{L of DPBS}$$

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 29 OF 64

Coat Chip-Arrays with ECM

Steps

1. Pour ECM solution into a disposable pipetting reservoir on ice to allow pipetting with multichannel pipette.
2. Using a multichannel P200 pipette, add 60 μ L of ECM solution to the bottom channel, pipetting through the bottom channel seeding port.
 - a. This volume is enough to fill the bottom channel as well as the bottom channel trough with ECM. The troughs should be full but not overflowing.
 - b. It is recommended to add ECM solution with 6 tips at a time using a multichannel pipette.
 - c. Double check that ECM flowed through all 12 Emulations for each Chip-Array. If there is no flow, this is likely due to an improper interface of pipette tip with the bottom channel seeding port, or there is misalignment of the pipette guide. Attempt to add ECM solution to the bottom channel using a single channel P200 pipette. If this still does not properly flow, then detach and reattach the pipette guide and attempt re-coating that specific Emulation. See Chip-Array Handling section for steps on pipette guide detachment and attachment.
3. Using a multichannel P200 pipette, add 60 μ L of ECM solution to the top channel, pipetting through the top channel seeding port.
 - a. This volume is enough to fill the top channel as well as the top channel trough with ECM. The troughs should be full but not overflowing.
 - b. It is recommended to add ECM solution with 6 tips at a time using a multichannel pipette.
 - c. Double check that ECM flowed through all 12 Emulations for each Chip-Array. If there is no flow, this is likely due to an improper interface of pipette tip with the bottom channel seeding port, or there is misalignment of the pipette guide. Attempt to add ECM solution to the top channel using a single channel P200 pipette. If this still does not properly flow, then detach and reattach the pipette guide and attempt re-coating that specific Emulation. See Chip-Array Handling section for steps on pipette guide detachment and attachment.
4. Repeat steps 2-3 for all Chip-Arrays.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 30 OF 64

5. Visually inspect the Emulations and wash any bubbles from the channel with the ECM solution.
6. To prevent evaporation during incubation, add PBS to the PBS humidity troughs between each Emulation (each PBS trough is filled with ~150 μ L).
7. Place pipette guide cover over the Emulations, ensuring that the cover is placed properly and flush. Keep this lid on whenever moving the consumable from the AVA, tissue culture incubator, and the biosafety cabinet. Failure to do so will compromise sterility.
8. Wrap a piece of Parafilm around edges of each Chip-Array and pipette guide cover to help maintain sterility when Chip-Arrays are incubating overnight.
9. Incubate all coated Chip-Arrays overnight at 4°C.
 - a. *Note:* Chip-Arrays can be stored at 4°C for up to 2 days if kept moist.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 31 OF 64

Day 0: Introduce Cells to Chip-Arrays

Overview

Goals

- Seed cells on top channel of Chip-Array

Required Materials

- 1X DPBS (at room temperature)
- Cell culture media (cell-specific)
- Serological pipettes
- Single channel and multichannel pipettes and filtered tips
- Aspirator and sterile aspirating tips
- 50 mL conical tubes
- 50 mL disposable pipetting reservoirs
- Diluted trypan blue counting solution
- Hemocytometer
- Nunc™ 96-Well Polystyrene Round Bottom Microwell Plates
- 70% ethanol
- Microscope
- Ice bucket and ice (if needed)

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 32 OF 64

Trypan Blue Cell Counting Solution

Reagent	Volume	Source	Cat. No.
Cell Culture Medium	40 μ L	-	-
Trypan Blue	5 μ L	Sigma	93595

- Maintain the Trypan Blue Cell Counting Solution at room temperature prior to use.
 - Prepare the Trypan Blue Cell Counting Solution fresh for each use.
-

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 33 OF 64

Prepare Chip-Arrays

Steps

1. Transfer all ECM-coated Chip-Arrays from 4°C to the AVA to incubate at 37°C for one hour.
 2. Fully aspirate the ECM solution from the top and bottom channel troughs, then from each channel by gently inserting aspirating pipette in top and bottom channel seeding ports. All Emulations should be dry after this step with no remaining ECM in the channels.
 3. Wash the bottom channel, then the top channel twice with 50 µL of warm cell culture media, slowly pipetting and aspirating the media coming out of the trough for each channel.
 - a. It is recommended to wash channels with 6 tips at a time using a multichannel pipette.
 - b. After the first wash, make sure to fully aspirate the troughs so that they are dry, while being careful not to aspirate media from the channels.
 - c. Hold the aspirator along the far edge of the trough and avoid the port when aspirating the top channel outflow.
 - d. After the second wash, leave the outflow volume of media in the top and bottom channel troughs to keep the Emulations hydrated until seeding.
 4. Thoroughly dry the area around the top channel seeding port with an aspirating pipette to prepare for cell seeding, being careful not to aspirate the channel.
 5. Cover the Emulations with the pipette guide cover and return Chip-Arrays to the AVA until the cells are ready for seeding.
-

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 34 OF 64

Prepare Cells

Overview

The preparation (thawing/harvesting) of cells is specific to the organ-model. It is essential to ensure the seeding density for your organ model is optimized for viable, functional cells and long-term culture. This protocol assumes the top channel is seeded first.

NOTE: This protocol assumes that top and bottom channel seeding will be done on Day 0 and Day 1, respectively.

Steps

1. Generate homogeneous cell solution in cell culture media
2. Measure the total suspension volume using a P1000 or 2 mL pipette.
3. Transfer 5 μ L of the homogeneous cell suspension to the Trypan Blue Cell Counting Solution, generating a 1:10 dilution.
4. Mix the trypan blue solution thoroughly and count the cells using a hemocytometer.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 35 OF 64

Cell Counting and Viability Assessment

- Count both viable and non-viable cells in each quadrant of the hemocytometer.

Live Cell Count, Dead Cell Count, Total Cell Count.

- Calculate the percentage viability of the cell solution.

Live Cells / Total Cells x 100 = % Viability

- Calculate the viable cell concentration. The dilution factor is 10 when prepared in the trypan blue solution above.

Live Cell Count x (1 x 10 x 10⁴) / 4 = Viable Cell Concentration (cells / mL)

- Calculate the viable cell yield.

Viable Cell Concentration x Cell Suspension Volume = Viable Cell Yield (cells)

Viable Cell Yield / Desired Density = Reconstitution Volume

Diluting Cells

After calculating the Viable Cell Yield, dilute cells with warm Medium to the required final cell density.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 36 OF 64

Seed Cells to the Top Channel

Overview

NOTE: This protocol assumes that top and bottom channel seeding will be done on Day 0 and Day 1, respectively. Prime Chip-Array will follow immediately after bottom channel cell incubation in this protocol. Users can modify the timeline of top and bottom channel seeding depending on the specific needs of the organ-model.

Work with one Chip-Array at a time per user. After seeding the first Chip-Array Emulation, use a microscope to assess the cell density within the channel. Adjust the density of the cell suspension as necessary for the rest of the chips.

Steps

1. Transfer the prepared Chip-Arrays from the AVA to the BSC.
 - a. When seeding, work with one Chip-Array at a time per user. Leave other Chip-Arrays in a 37°C incubator or the AVA.
2. Carefully aspirate media from the top channel troughs for all Emulations, ensuring not to touch the trough port linked to the top channel.
 - a. **NOTE:** Having a completely dry top channel trough before seeding is essential in ensuring uniformity and even distribution of cells in the top channel. Having any liquid in the top channel trough before seeding may encourage movement of cells due to hydrostatic pressure.
3. Designate the first Emulation on a Chip-Array to test the seeding density.
4. Very gently agitate the cell suspension before seeding to ensure a homogeneous cell suspension.
5. Quickly and steadily pipette 20 µL of cell solution with a single channel P20 pipette in the top channel through the top channel seeding port, aspirating the outflow from the top channel trough immediately without touching the trough port.
 - a. When seeding, it is recommended to leave the P20 pipette tip in the seeding port until the outflow volume in the trough is aspirated, then the user may remove the pipette tip.
6. Cover the Chip-Array with the pipette guide cover and visually inspect the seeded Emulation under a lab microscope to check if the density and distribution look appropriate (see Figure 13 below). Cells should not be overcrowded and

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 37 OF 64

should have an even layer with ~one cell radius between individual cells.
Troubleshoot by readjusting density if not.

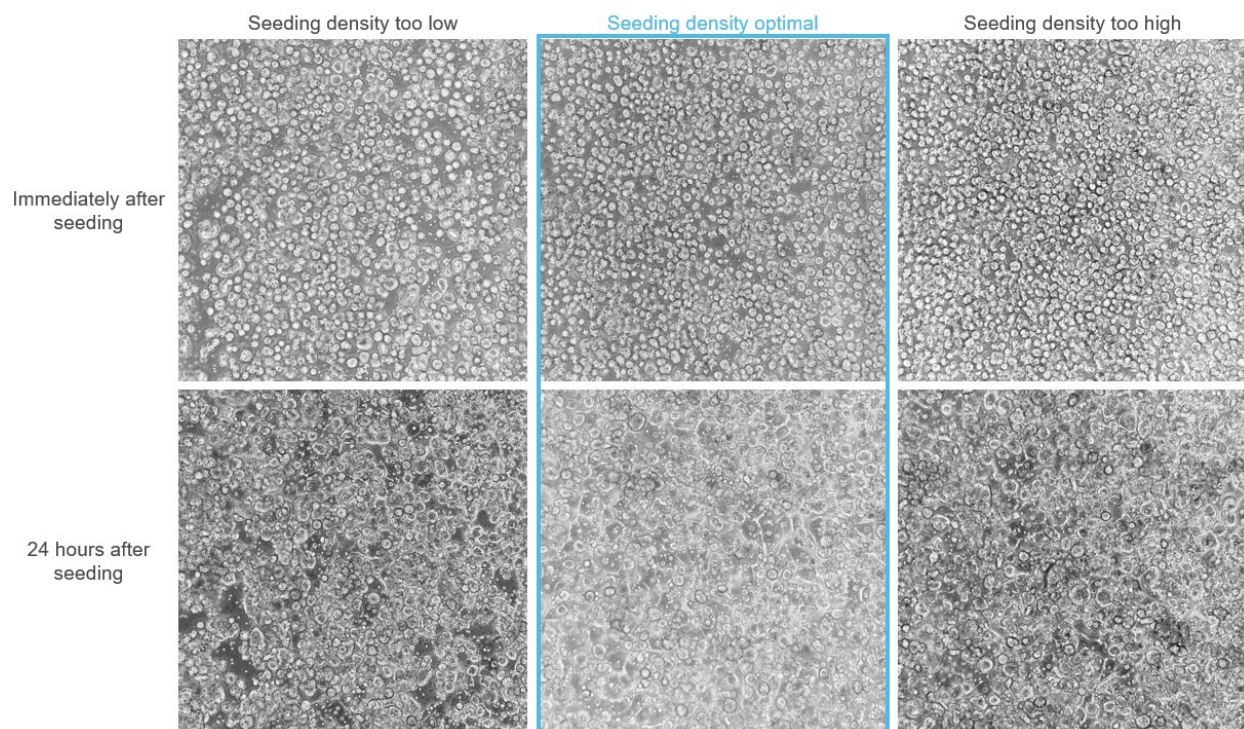


Figure 13. Example seeding density reference chart (Hepatocytes, Liver Quad-Culture)

- If cell distribution is not uniform due to gaps or gradients present in the top channel, vacuum aspirate the top channel of the specific Emulation, then flush the top channel twice with 50 μ L Media while aspirating the outflow.
- Re-seed the Emulation again with 20 μ L cell suspension with a single channel P20 pipette and double check cell distribution.
- If the seeding density is too dense or not optimal, wash the test Emulation according to Step 6a, adjust the cell density accordingly, and re-seed the test Emulation until the density in the channel is correct.
- If excessive fluid flow is observed during cell settling (referred to as “cell racing”), an extra step is required to ensure that cells settle on the top channel membrane evenly. Add 500 μ L of Media to the top outlet reservoir of the Emulation experiencing cell racing. Then proceed with the washing and re-seeding steps as described in Steps 6a and 6b, respectively.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 38 OF 64

7. After confirming the correct cell density, agitate the cell suspension to ensure a homogenous cell suspension, then transfer 400 μ L of cell solution to 6 wells of a 96 well Ultra-Low Attachment (ULA) plate.
8. Resuspend the cell suspension in the wells using a multichannel P200 pipette.
 - a. If you are working with cells that readily settle to the bottom of the wells, it is recommended to resuspend the cell suspension in these wells before seeding each Chip-Array.
9. On the same Chip-Array, seed Emulations 2-6, then Emulations 7-12 with a multichannel P20 pipette, pipetting quickly and steadily 20 μ L of cell solution through the top channel seeding port. Allow cells to settle for a minute before checking for appropriate distribution using a lab microscope.
 - a. *NOTE:* When seeding several Emulations at a time, users must be careful with the amount of force applied to the multichannel P20 tips. Inserting the tips is enough to create a seal with the chip gasket to ensure proper flow. Additional excessive force when pipetting multiple Emulations at a time (i.e. using your offhand to apply force to the multichannel pipette) can cause tips to become stuck in the pipette guide.
 - b. Always make sure tips are perpendicular when inserted into the pipette guide; failure to do so can cause tips to become stuck.
 - c. It is recommended to seed no more than 6 Chip-Array Emulations at a time. Seeding more than 6 Emulations simultaneously can potentially lead to variability and pipetting issues.
10. After confirming the correct cell density and seeding the first Chip-Array, seed all remaining Chip-Arrays using a multichannel P20 pipette, working in groups of 6 Emulations at a time.
 - a. *Note:* Minimize the amount of time the cells are outside the incubator by seeding one Chip-Array at a time per user and then immediately placing each seeded Array into the AVA at 37°C or a 37°C incubator.
11. Capture representative images on a lab microscope (3-4 Emulations per Chip-Array).
 - a. If any Emulations need to be reseeded, follow the protocol outlined in Step 6
12. If needed, top off the bottom channel trough with warm Media to ensure the bottom channel does not dry out during incubation.
13. Add any additional PBS to the PBS troughs if needed.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 39 OF 64

14. Incubate Chip-Arrays in the AVA at 37°C for 4 hours. See Figure 14. for examples of attachment.
15. After incubation, **gently** wash top channel once with 50 μ L of cell culture media to clear away debris and non-attached cells. **Gently** wash top channel again and leave cell culture media in top channel trough to prevent evaporation.
16. If seeding bottom channel cells immediately, proceed to next section. If not, incubate Chip-Arrays overnight in 37°C incubator.

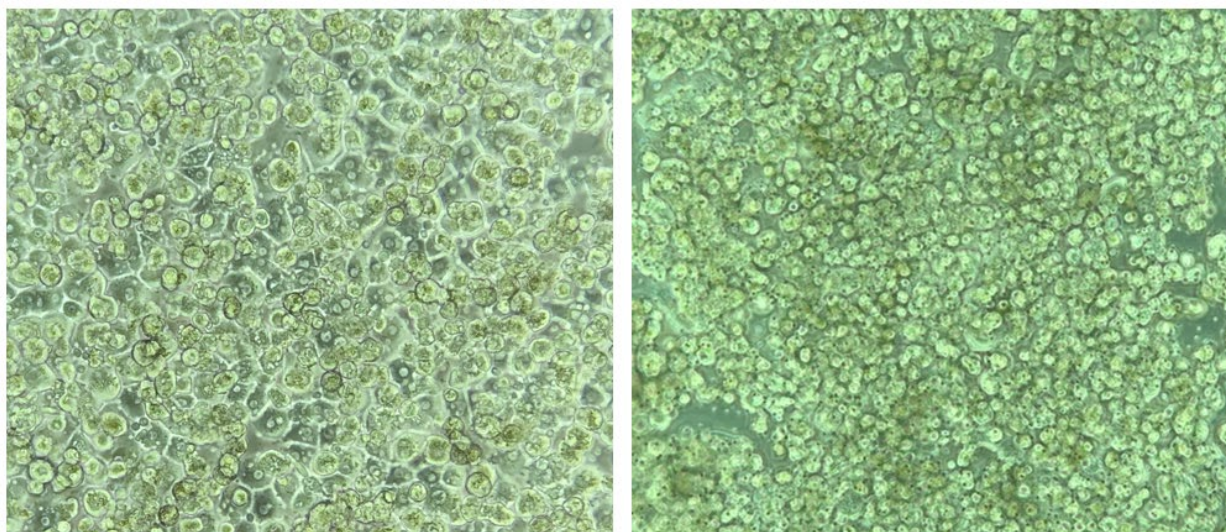


Figure 14. Example of appropriate (left) and poor (right) cell attachment (Hepatocytes, Liver Quad-Culture)

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 40 OF 64

Day 1: Chip-Array Prime, Regulate, and Flow

Overview

Goals

- Wash Chip-Arrays
 - Seed cells on bottom channel of Chip-Array
 - De-gas and equilibrate media
 - Prime Chip-Arrays
 - Regulate Chip-Arrays and begin flow
-

Materials

- Trypan blue counting solution
- Hemocytometer
- 50 mL conical tube
- Serological pipettes
- Single channel and multichannel pipettes and filtered tips
- Aspirator and sterile tips
- 50 mL disposable pipetting reservoir
- Nunc™ 96-Well Polystyrene Round Bottom Microwell Plates
- Ice bucket and ice
- 70% ethanol
- Microscope
- Installed and qualified AVA instrument
- Steriflip® filtration unit: PVDF filter 0.45 µm (sterile)

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 41 OF 64

Wash Chip-Arrays

Steps

1. Prepare Medium and warm to 37°C in bead bath. Aliquot Medium to keep cool at 4°C.
2. Image representative Emulations to capture cell morphology 24h after seeding.
3. **Gently** wash bottom channel by pipetting twice with 50 µL of warm Medium through the bottom channel seeding port, aspirating the media coming out of the trough and being careful not to aspirate media from the channel.
4. **Gently** wash top channel by pipetting twice with 50 µL of warm Media through the top channel seeding port, aspirating the media coming out of the trough and being careful not to aspirate media from the channel.
5. Aspirate PBS from PBS troughs and refill each PBS trough to exactly 100 µL. This is enough volume to partially fill the troughs while also not spilling when inverting the Chip-Array.
6. Use previous techniques for multichannel P200 pipette and ensure port area and bottom channel trough is dry for seeding. Return the Chip-Arrays to the incubator.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 42 OF 64

Seed Cells to the Bottom Channel

Overview

NOTE: This protocol assumes that top and bottom channel seeding will be done on Day 0 and Day 1, respectively. Prime Chip-Array will follow immediately after bottom channel cell incubation in this protocol. Users can modify the timeline of top and bottom channel seeding depending on the needs of the organ-model.

Repeat procedure from “Prepare Cells” and “Cell Counting and Viability Assessment” for bottom channel cells.

Steps

1. Transfer the prepared Chip-Arrays from the AVA to the BSC.
2. Carefully aspirate media from the bottom channel troughs for all Emulations, ensuring not to touch the trough port linked to the bottom channel.
 - a. **NOTE:** Having a completely dry bottom channel trough before seeding is essential in ensuring uniformity and even distribution of cells in the bottom channel. Having any liquid before seeding may encourage movement of cells due to hydrostatic pressure.
3. Designate the first Emulation on a Chip-Array to test the seeding density.
4. Very gently agitate the cell suspension before seeding to ensure a homogeneous cell suspension.
5. Quickly and steadily pipette 10 μ L of cell solution with a single channel P20 pipette in the bottom channel through the bottom channel seeding port, aspirating the outflow from the bottom channel trough immediately without touching the trough port.
 - a. When seeding, it is recommended to leave the P20 pipette tip in the seeding port until the outflow volume in the trough is aspirated, then the user may remove the pipette tip.
6. Cover the Chip-Array and quickly visually inspect under a lab microscope if the density and distribution look appropriate (see Figure 15). Troubleshoot by readjusting density if not.
 - a. If distribution isn't uniform due to gaps or gradients present in the bottom channel, flush the bottom channel with 50 μ L cell culture media twice,

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 43 OF 64

while aspirating the outflow. Do not vacuum aspirate the bottom channel, as this may remove media from the top channel.

- b. Re-seed the Emulation again with 10 μ L cell suspension with a single channel pipette.
- c. If the seeding density is not optimal, wash the test Emulation according to Step 6a. above, adjust the cell density accordingly, and re-seed the test Emulation until the density in the channel is correct.



Figure 15. Example of optimal cell seeding density in bottom channel (NPCs, Liver Quad-Culture)

7. After confirming the correct cell density, agitate the cell suspension to ensure a homogenous cell suspension, then transfer 150 μ L to 6 wells in ULA plates.
 - a. Place ULA plate on ice to keep the cells cool, if needed
8. Resuspend the cell suspension in the wells using a multichannel P200 pipette.
 - a. Take care while resuspending as such a small volume is prone to bubbles that reduce the total volume available.

- b. If left static for too long, the cells will settle to the bottom of the wells. It is recommended to resuspend the cell suspension in these wells with a multi-channel pipette before seeding each Chip-Array.
9. On the same Chip-Array, seed Emulations 2-6, then Emulations 7-12 with a multichannel P20 pipette, pipetting quickly and steadily 10 μ L of cell solution through the bottom channel seeding port. Invert Array as swiftly as possible.
 - a. *NOTE:* When seeding several Emulations at a time, users must be careful with the amount of force applied to the multichannel P20 tips. Inserting the tips is enough to create a seal with the chip gasket to ensure proper flow. Additional excessive force when pipetting 6 Emulations at a time (i.e. using your offhand to apply force to the multichannel pipette) can cause tips to become stuck in the pipette guide.
 - b. Always make sure tips are perpendicular when inserted into the pipette guide; failure to do so can cause tips to become stuck.
 - c. Be very careful when inverting Chip-Arrays, making sure that hands are gripping both the pipette guide cover and the reservoir lid. Failure to do so can cause either of the lids to fall off, potentially compromising sterility.
10. Seed all remaining Chip-Arrays using a multichannel P20 pipette working in groups of 6 Emulations at a time, imaging representatives immediately and inverting the Chip-Array as swiftly as possible.
 - a. If any Emulations need to be reseeded, flush the bottom channel with 50 μ L cell culture media twice, while aspirating the outflow. Re-seed the Emulation again with 10 μ L cell suspension with a single channel pipette.
 - b. *Note:* Minimize the amount of time the cells are outside the incubator by seeding one Chip-Array at a time per user and then immediately placing each seeded, inverted Array into an incubator at 37°C.
11. Incubate inverted Chip-Arrays in an incubator at 37°C for 4 hours.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 45 OF 64

Gas Equilibration of Media

CAUTION: The media equilibration step is imperative to successfully culture Organ-Chips. Omitting this step will cause bubbles to form in the Chip-Array, which will in turn negatively impact cell viability and can result in irregular flow.

Before You Begin

- Once the culture media for both chip channels have been warmed for at least 1h, the gas equilibration step can be performed.
- Ensure the medium is outside of a warmed environment (such as an incubator or bath) for no longer than 10 minutes, as gas equilibrium can become compromised when medium is allowed to cool.
- If the vacuum pump is not located close to the water bath (e.g., inside the BSC), it is recommended to place some clean water warmed at 37°C inside the BSC to minimize cooling during the media equilibration step.

Steps

1. Prepare and warm cell culture media
 - a. Once the culture media has been warmed for at least 1 hour, the gas equilibration step can be performed.
2. Place at least 2 mL of Medium for each Emulation in a 50 mL conical tube.
3. Perform step 2 for medium to be used in the experiment.
4. Warm all conical tubes of media at 37°C in water bath or bead bath for at least 1 hour.
5. Immediately connect the 50 mL tubes containing each warmed medium to a Steriflip unit using the following steps:
 - a. Attach each conical tube containing warmed media to a Steriflip unit.
 - b. With the unit “right-side up” (medium in the bottom conical tube), apply vacuum and wait for 10 seconds.
 - c. Invert the Steriflip-connected tubes, and check that the medium begins to pass from the top conical tube to the bottom tube.
 - i. *Note:* The vacuum must operate at a minimum of -70 kPa. At this pressure, it should take about 2 seconds for 10 mL of media to flow

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 46 OF 64

through the filter. If it takes longer, stop and refer to “Media taking too long to pass through Steriflip” in the troubleshooting section.

- d. Leave the filtered medium under vacuum for at least five minutes.
6. Remove the vacuum tubing from the Steriflip unit.
7. Separate the conical tubes containing media from the Steriflip unit and immediately place them into the incubator with the caps loose to maintain the dissolved gas equilibrium state and allow bubbles to escape.
 - a. *Note:* Minimize the time media is outside of the incubator when the Chip-Arrays are being prepared to maintain the correct temperature. This is critical to ensure Chip-Array success.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 47 OF 64

Prime Chip-Arrays

CAUTION: Ensuring successful Chip-Array priming is best practice for successful liquid-liquid connection between the Emulation and the Chip-Array. Failure to do so can cause bubbles to form in the Chip-Array, which will in turn negatively impact cell viability and result in irregular flow.

Steps

For the following steps, handle Chip-Arrays one at a time, setting each Chip-Array aside in the incubator after media is added to reservoirs before Prime:

1. After incubating bottom channel cells for 4 hours, return Chip-Arrays to the “right side up” by flipping consumable over.
 - a. Be very careful when flipping Chip-Arrays, making sure that hands are gripping both the pipette guide cover and the reservoir lid. Failure to do so will cause either of the lids to fall off, potentially compromising sterility.
2. Image several representative Emulations to ensure proper attachment is observed. See Figure 16. for examples of attachment in the bottom channel.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 48 OF 64

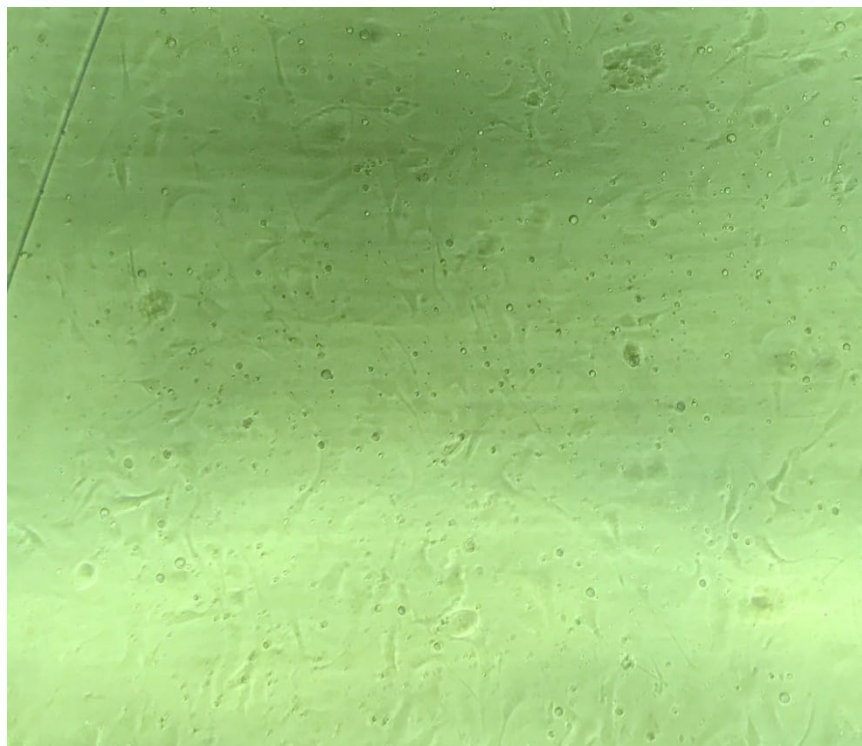


Figure 16. Example of optimal cell attachment in bottom channel (NPCs, Liver Quad-Culture)

3. Gently wash both top channel and bottom channel with 50 μ L of their respective warmed and equilibrated media, leaving a small amount of media in the pipette tips after dispensing. Allow media to fill the pipette guide troughs – do not aspirate while dispensing media.
 - a. Do not fully dispense media from the pipette tips. Leaving a small amount of media in the pipette tip rather than dispensing the full volume helps ensure large bubbles are not trapped in the seeding ports.
4. Aspirate media from the pipette guide troughs and humidity troughs (PBS filled).
 - b. Place aspirating pipette tip in the corner of the troughs and avoid contact or proximity to the port leading to the emulations.
 - c. Allow media to be fully aspirated from the troughs before moving to the next trough for aspiration. This may take over one second to achieve.
5. Place the pipette guide cover back over emulations.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 49 OF 64

6. Remove the tape tabs from the sides of the consumable and peel the plastic protector from the reservoir lid.
7. Fill the top channel inlet reservoirs by adding 1000 μ L of warmed and equilibrated media directly over the top channel inlet vias using a multichannel P1000 pipette.
 - d. Reservoir vias are located in the bottom left corner of each reservoir compartment. The recommended technique is to hug the bottom left corner of the inlet reservoirs with the pipette tips, then allow liquid to dispense down the wall onto the via – this minimizes the risk of trapping air pockets in the vias.
 - e. See Figure 17 for locations of reservoir vias.
8. Fill the top channel outlet reservoirs by adding 500 μ L of warm and equilibrated media directly over the top channel outlet vias using a multichannel P1000 pipette.
9. Fill the bottom channel inlet reservoirs by adding 1000 μ L of warmed and equilibrated media directly over the bottom channel inlet vias using a multichannel P1000 pipette.
 - f. Reservoir vias are located in the bottom left corner of each reservoir compartment. The recommended technique is to hug the bottom left corner of the inlet reservoirs with the pipette tips, then allow liquid to dispense down wall onto the via – this minimizes the risk of trapping air pockets in the vias.
10. Fill the bottom channel outlet reservoirs by adding 500 μ L of warmed and equilibrated media directly over the bottom channel outlet vias using a multichannel P1000 pipette.
11. The final volume of media in the reservoirs at this step:
 - g. All inlet reservoirs should be filled with 1000 μ L of media.
 - h. All outlet reservoirs should be filled with 500 μ L of media.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 50 OF 64

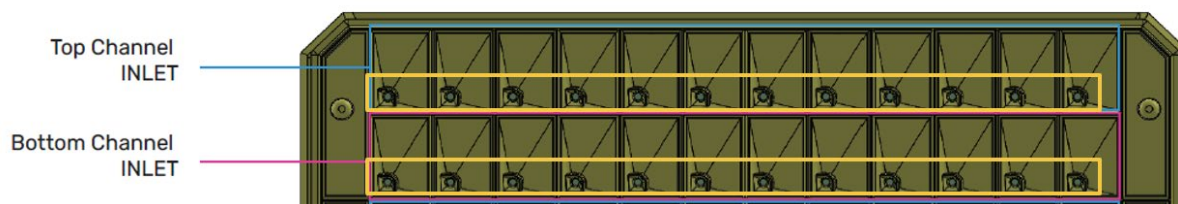


Figure 17. Location of reservoir vias.

12. Place the reservoir lid on top of the reservoirs. Minimize any motion that may splash media onto the reservoir lid.
13. Remove pipette guides by pushing on the clip then lifting upwards. Once the clip is at a 45-degree angle, lightly tug on the clip to remove it. Be careful not to drag your hand over media in the gasket ports.
 - a. Refer to “Attachment and Detachment of Pipette Guides and Flow Guides onto Chip-Array” section for more guidance on pipette guide removal.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 51 OF 64

14. Aspirate any excess media from the viewing windows and on the gasket. Do not aspirate within the gasket ports. See Figure 18 for proper aspiration of excess media around the gasket ports.

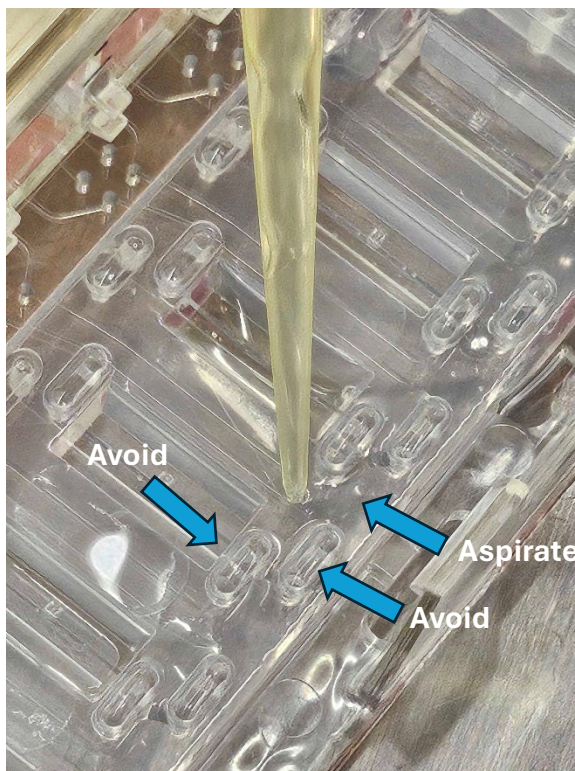


Figure 18. Proper aspiration of excess media around Emulation gasket ports.

15. Place the pipette guide cover back onto the Chip-Array.
 - a. Do this one by one for all Chip-Arrays, setting each Chip-Array aside in the incubator when media has been added to the reservoirs.
 - b. Once all Chip-Arrays are ready for prime, bring them out of the incubator and to the AVA. Be careful not to tip or tilt the Chip-Arrays as this will cause media to splash onto the reservoir lids.
16. Eject the trays using the AVA tray button or using the AVA software.
17. Place the Chip-Arrays on the trays, making sure that the Emulations are facing inwards towards the AVA instrument. Double check that each Chip-Array is slotted flush against the pins on each tray.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 52 OF 64

- a. **Additionally, ensure that all reservoir lids are flat and secure on the Chip-Arrays before inserting the trays.** This can be checked by pushing forward on the corners of the lid facing the user as the Chip-Arrays sit on the AVA trays. Afterwards, press the lid flat. These steps will ensure that the lids are flush with the Chip-Arrays. Failure to perform this check can lead to flow issues. **Perform this check every time the Chip-Arrays are inserted into the AVA.**
18. Insert the trays using the AVA tray button or using the AVA software.
19. Click the “Flow Operations” dropdown menu and select Prime. In the box below the Prime button, the percent progress will appear.
 - a. Prime Cycle parameters (Inlets 30 kPa, Outlets 0 kPa for 45 seconds).
 - b. The Prime Cycle is performed to fill the microfluidic channels on the inlet and outlet sides of the Chip-Array Emulations with media from the Chip-Array inlet reservoirs.
20. After Prime Cycle has completed, eject the trays using the AVA tray button or using the AVA software.
21. Return to BSC and double check that priming was successful by observing media filling the gasket port cavities. See Figure 19 for an example of successful priming.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 53 OF 64



Figure 19. Successful priming indicated by media visible in the gasket port cavities.

22. “Flood” the inlet gasket ports by adding large droplets of media over the inlet gasket ports. This should create large pools of media over the inlet gasket ports. Any bubbles in the gasket ports post-Prime should now rise to the surface of these large pools of media.
23. Aspirate any excess bubbles on top of the large droplets of media around the gasket port cavities, ensuring that domed droplets still remain above the inlet gasket ports. Aspirate any excess media in the viewing windows of the Chip-Array. Do not aspirate within the gasket ports.
 - a. The outlet port cavities may have a large droplet escaping the cavity. Typically, the media will be contained within the cavity and will either be level with the top of the gasket or protrude slightly above the top of the gasket.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 54 OF 64

- b. Be careful to avoid aspirating media out of the top and bottom channels. Do not aspirate media from within the gasket ports.
- 24. Attach the flow guides to the Chip-Arrays, **but do not snap them closed. Leave the flow guides open at ~45° angle.**
 - a. When attaching flow guides, a light “click” sound should be heard by the user when the guide is connected to the reservoir hinge.
 - b. Once flow guides are attached, leave them open at an ~45° angle and do not push them closed downwards.
 - c. Refer to “Attachment and Detachment of Pipette Guides and Flow Guides onto Chip-Array” section for more guidance on flow guide attachment and snapping into place.
- 25. Tilt Chip-Arrays forward at a 45° angle for 10 seconds, or until multiple large droplets form at the outlet gasket ports. Place the Chip-Array flat on the bench for 2 seconds, and then gently snap the flow guides in place one at a time, using two fingers to hold onto the flow guide tab while pressing downward. This process ensures consistent priming of the emulations.
 - a. A loud click should be heard as the flow guides are snapped closed.
 - b. Once the flow guides are connected to the Emulations, this creates a seal that prevents exposure of the Emulations to the outside environment, helping to ensure sterility. After flow guides are attached, there is no need to use the pipette guide cover anymore. These can be set aside.
 - c. Double check that all flow guides are correctly snapped into place by pressing onto the flow guides.
 - d. Double check the reservoir lid. If excess media is observed on the lid gasket, vacuum aspirate this excess media.
- 26. Move all Chip-Arrays to the AVA.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 55 OF 64

Regulate Chip-Arrays

Steps

1. Eject the trays using the AVA tray button or using the AVA software.
2. Place the Chip-Arrays on the trays, making sure that the Emulations are facing inwards towards the AVA instrument. Double check that each Chip-Array is slotted flush against the pins on each tray.
3. Insert the trays using the AVA tray button or using the AVA software.
4. Select the “Flow Operations” dropdown menu, select Regulate, select which manifolds should be lowered based on Chip-Array position in the AVA, and set flow rate parameters that the system will default to after Regulate is complete. Top and bottom channel flow rates should be set to the recommended flow rate for your organ model.
5. Manifolds will lower and Regulate will begin. In the box below the Regulate button, the percent progress will appear.
 - a. Regulate Cycle parameters:
 - i. Total Duration: ~4 hours

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 56 OF 64

Day 2: Chip-Array Regulate

Overview

Goals

- Perform 2nd Regulate on Chip-Arrays
- Maintain Chip-Arrays in AVA

Required Materials

- Chip-Arrays
- Cell culture media

Regulate Chip-Arrays

Steps

1. Prepare and warm cell culture media. Once media has been warmed for at least 1 hour, the gas equilibration step can be performed according to the “Gas Equilibration of Media” section above.
2. Pause flow on the AVA and transfer two Chip-Arrays at a time from the AVA to the BSC.
3. Aspirate all inlet and outlet reservoirs.
4. Replenish inlet reservoirs with 1250 μ L of respective, equilibrated media and outlet reservoirs with 500 μ L of respective, equilibrated media.
5. Eject the trays using the AVA tray button or using the AVA software.
6. Place the Chip-Arrays on the trays, making sure that the Emulations are facing inwards towards the AVA instrument. Double check that each Chip-Array is slotted flush against the pins on each tray.
7. Insert the trays using the AVA tray button or using the AVA software.
8. Select the “Flow Operations” dropdown menu, select Regulate, select which manifolds should be lowered based on Chip-Array position in the AVA, and set flow rate parameters that the system will default to after Regulate is complete. Top and bottom channel flow rates should be set to the recommended flow rate for your organ model.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 57 OF 64

9. Manifolds will lower and Regulate will begin. In the box below the Regulate button, the percent progress will appear.
 - a. Regulate Cycle parameters:
 - i. Total Duration: ~4 hours
10. It is recommended after the 2nd Regulate is performed to remove all Chip-Arrays, aspirate the outlets, and top off the media in the inlets up to 1250 μL to facilitate flow over the next three days (if flow rate is set to 15 $\mu\text{L}/\text{h}$).

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 58 OF 64

Day 5+: Chip-Array Maintenance and Sampling

Overview

Goals

- Maintain Chip-Arrays in AVA
- Cell culture inspection
- Collect samples for analysis

Required Materials

- Chip-Arrays
- Cell culture media

CAUTION: Failure to replenish media on time will cause air to be introduced into the Chip-Array, which ultimately results in cell death. Filling inlet reservoirs up to 1250 μL will ensure up to 72 hours of flow, if flow rate is set to 15 $\mu\text{L}/\text{h}$. Do not fill the inlet reservoirs past 1300 μL of the total volume during replenishment.

Changing Media

Steps

1. Media should be refilled every 2 days to ensure the inlet reservoirs do not deplete of media. At 15 $\mu\text{L}/\text{h}$, the maximum time between media replenishments should be 72 hours.
2. If compound dosing is part of your experiment design (see Dosing section), make sure the proper dosing media is added to the correct Chip-Array Emulations.
 - a. Warm cell culture media for 1 hour and dose with compound if necessary.
3. Pause flow on the and transfer two Chip-Arrays at a time from the AVA to the BSC.
4. Remove Chip-Array lid and collect effluent from the Chip-Array inlet and outlet reservoirs, taking care not to disturb the reservoir vias.
5. Any media not collected for analysis should be aspirated by aspirating from the top right corner of the reservoirs farthest from the reservoir via while the

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 59 OF 64

consumable lies flat. This ensures that a thin liquid film still covers the reservoir Vias so that no air is introduced into the fluidic channels of the Chip-Array.

- a. When aspirating reservoirs, it is recommended to use a multichannel aspirator to streamline the aspirating workflow. This will help to prevent cross-contamination of dosing media across different conditions. Users should also prevent cross-contamination of top and bottom channel media when aspirating.
6. Refill the Chip-Array inlet reservoirs with 1250 μ L of respective top and bottom channel warmed media.
7. Place the Chip-Array reservoir lids on and transfer the Chip-Arrays back to the AVA (two at a time) to resume flow.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 60 OF 64

Automated Imaging

Steps

Automated imaging protocols should be assigned for every day required. The AVA software allows users to exclude Emulations if an Emulation is deemed unsuitable for assigning to a condition.

1. Users should not eject the trays or remove the Chip-Arrays when an automated imaging protocol is running. For this reason, it is recommended that users schedule Chip-Arrays to be imaged during specified times when the duration of the imaging does not interfere with lab activities.
2. It is recommended that at least 4 fields of view (FOVs) be captured for every Emulation.
3. After each automated imaging session, these image sets can be viewed in the Image Reviewer tab.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 61 OF 64

Dosing

Steps

1. Assign control and dosing conditions to Chip-Array Emulations.
2. Aspirate media from the Chip-Array reservoir inlets and outlets (handle Chip-Arrays two at a time to minimize time outside of AVA or BSC).
3. Fill the reservoir inlets with their corresponding dosing media based on the condition assigned to the specific Emulation.

Sampling

Steps

1. Effluent can be collected from the Chip-Array outlet reservoirs and transferred to a collection plate for analysis.
 - a. It is recommended to handle consumables two at a time when sampling, returning consumables to the AVA with flow set to "0" immediately after taking samples.
2. Remove Chip-Array lid and collect effluent from the Chip-Array inlet and outlet reservoirs, taking care not to disturb the reservoir vias.
 - a. *Note:* It is recommended to freeze the effluent at -80°C if effluent will not be analyzed immediately. It is also recommended to reduce the number of freeze-thaw cycles to prevent degradation of proteins or cytokines in the effluent, such as albumin.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 62 OF 64

Troubleshooting

Bubbles and Flow Issues

If a bubble is present in the channel during ECM coating, washing, or cell seeding, it is recommended to wash the channel with the appropriate solution until all bubbles have been removed. If bubbles persist, it may be helpful to aspirate the channel dry and slowly re-introduce solution.

After flow has been initiated on the AVA, it is possible that users may run into issues related to bubble formation and blockage in the fluidic elements of the consumables. If these are severe enough, they may cause lower than expected flow or no flow at all. This can usually be detected by eye when assessing volume in the top or bottom channel outlets. If this occurs, this can potentially be resolved via numerous methods based on the issue:

1. If a bubble is observed visually in the top and/or bottom channel, this bubble potentially can be removed by:
 - a. Flushing the Chip-Array at 1,000 $\mu\text{L}/\text{h}$ for 5 minutes. This should push the bubble out of the channel.
 - b. If Step 1a does not work and bubble is still present, then detach the flow guide for the specific Emulation, reattach the pipette guide, and manually flush the bubble out with a pipette.
 - i. Afterwards, it is highly recommended to detach the pipette guide, re-prime the Chip-Array, reattach the flow guide, and re-run Regulate on the Chip-Array.
 - ii. Due to this manual intervention, b. should be considered a last resort if a. does not work.
2. If a bubble is not observed visually in the top and/or bottom channel, this indicates the bubble is likely in the fluidic elements of the consumable. This bubble can potentially be removed by:
 - a. Flushing the Chip-Array at 1,000 $\mu\text{L}/\text{h}$ for 5 minutes. This should push the bubble out of the channel.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 63 OF 64

- b. If the flow issue persists afterwards, re-run Regulate on the Chip-Array.
There is no need to manually detach the flow guide for manual flushing with a pipette.
3. If media appears to be collecting in the outlet reservoirs during sampling, this could indicate the Chip-Array reservoir lid has media splashed on it and media is wicking out of the reservoirs. To mitigate, remove the reservoir lid and wipe clean with a cleanroom wipe sprayed with ethanol. Do not fill reservoirs beyond 1mL during media replenishment. Handle the Chips-Arrays in such a way so as to minimize inadvertent contact of media with the reservoir lid (e.g. tilting the Chip-Array during transport). Minimize time consumables are outside the AVA.

Media Taking Too Long to Pass Through Steriflip

If media takes too long to pass through Steriflip unit when equilibrating the dissolved gas in the media before Prime, it is likely that vacuum pressure is not reaching -70kPa. It is recommended to find an alternate vacuum source with the appropriate pressure.

Priming Issues

If Chip-Arrays do not properly prime, ensure that media are covering the reservoir Vias and run the Prime Cycle again on the AVA.

Environmental Controls

The AVA is set to maintain temperature and %CO₂ concentration with tolerances around these set points. If the AVA instrument is outside of these tolerances, users will be notified via an alert on the external workstation UI.

If these issues persist, contact Emulate Support to troubleshoot and diagnose root cause.

TITLE Chip-Array Basic Research Kit Protocol	DOCUMENT EP-244	REVISION B
	DATE 24Aug2026	PAGE 64 OF 64

Automated Imaging

The AVA software allows users to schedule “Automated Imaging” sessions, where the AVA microscope will image specific FOVs on specific Emulations designated by the user at a certain time/date.

If the AVA fails to capture these images, attempt to open a “Live Imaging” session and manually capture images to determine if this is a microscope failure.

If “Live Imaging” works, then this means the microscope is functional and the failed “Automated Imaging” session is likely to be related to the scheduling software. Refer to the AVA software manual or contact Emulate Support for further troubleshooting.